



Rapid Testing. Reliable Results.

Legionella Sampling Packing Instructions

Small Box (White Kraft 8W x 8D x 10.5H)

1 Sealable Insulated Bag

Option 1: Maximum of 18–250 mL Bottles

Option 2: Maximum of 4–1L Bottles

Option 1



Option 2



Open Box



Closed Box



Medium Box (Plain Kraft 16W x 16D x 9H)

1 Sealable Insulated Bag

Option 1: Maximum of 72–250 mL Bottles (18 per bag)

Option 2: Maximum of 16–1L Bottles (4 per bag)

Option 1



Option 2



Open Box



Closed Box



Large Bulk à la carte (28W x 18D x 18H)

of Small Boxes (White Kraft) – flat packed

of Medium Kit Boxes (Plain Kraft) – flat packed

of Sealable Insulated Bags – flat packed

of 250 mL Bottles

of 1L Bottles

Packing Instructions (Please Follow Carefully)

1. Insert sealable insulated bag into box
2. Load bag with bottles (250 mL and/or 1 L)
3. Remove cellophane and seal insulated bag
4. If headspace exists, fill the gap with adequate packing materials for secure shipment
5. Print off a copy of Chain of Custody/Sample Submission Form and place in package
6. Securely tape box closed for shipment
7. Attach shipping label (UPS/FedEx)
8. Transport box to carrier location (UPS/FedEx)